

## Gateway<sup>®</sup> BP Clonase<sup>™</sup> II Enzyme Mix

Cat. No. 11789-020

Size: 20 reactions


Cat. No. 11789-100

Size: 100 reactions

Store at -20°C (non-frost-free freezer)

### Gateway<sup>®</sup> Technology

The Gateway<sup>®</sup> Technology is a universal cloning method that takes advantage of the site-specific recombination properties of bacteriophage lambda (1) to provide a rapid and highly efficient way to move DNA sequences into multiple vector systems. The Gateway<sup>®</sup> Technology is schematically represented below.

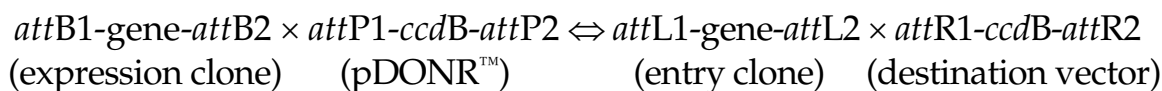

The *attB* × *attP* reaction is mediated by Gateway<sup>®</sup> BP Clonase<sup>™</sup> II enzyme mix; the *attL* × *attR* reaction is mediated by Gateway<sup>®</sup> LR Clonase<sup>™</sup> II enzyme mix. *ccdB* is the F plasmid-encoded gene that inhibits growth of *E. coli* (2,3) and “gene” represents any DNA segment of interest (e.g. PCR product, cDNA, genomic DNA).

### Description

Gateway<sup>®</sup> BP Clonase<sup>™</sup> II enzyme mix is a proprietary enzyme and buffer formulation containing the bacteriophage lambda recombination protein Integrase (Int), the *E. coli*-encoded protein Integration Host Factor (IHF) (1), and reaction buffer provided in a single mix for convenient reaction set up. Gateway<sup>®</sup> BP Clonase<sup>™</sup> II enzyme mix catalyzes *in vitro* recombination between an *attB*-PCR product (or *attB*-containing expression clone) and an *attP*-containing donor vector to generate an *attL*-containing entry clone. Store Gateway<sup>®</sup> BP Clonase<sup>™</sup> II enzyme mix at -20°C (non-frost-free freezer) for up to 6 months. For long-term storage, store at -80°C.

### Components Supplied

Components Supplied	20 rxns	100 rxns
Gateway <sup>®</sup> BP Clonase <sup>™</sup> II Enzyme Mix	40 µl	200 µl
Proteinase K Solution (2 µg/µl)	40 µl	200 µl
30% PEG 8000/30 mM MgCl <sub>2</sub> Solution	1 ml	5 ml
pEXP7-tet Positive Control (50 ng/µl)	20 µl	20 µl

### Quality Control

BP Clonase<sup>™</sup> II enzyme mix is functionally tested in a 1 hour recombination reaction followed by a transformation assay.

Part No. 11789.II.pps

Rev. Date: 22 July 2005

## General Recommendations and Guidelines

- pEXP7-tet is provided for use as a positive control in the BP reaction and contains an approximately 1.4 kb fragment consisting of the tetracycline resistance gene and its promoter flanked by *attB* sites.
- For *attB*-containing expression clones, we recommend using plasmid DNA purified with the PureLink™ HQ Mini Plasmid Purification Kit (Catalog no. K2100-01). Mini-prep (alkaline lysis) DNA preparations are adequate for Gateway® cloning reactions; however, in general, such DNA cannot be quantitated by UV absorbance due to contaminating RNA and nucleotides.
- You may use *attB*-PCR products in the BP reaction without purification. To achieve a higher percentage of desired clones, use PEG/MgCl<sub>2</sub> precipitation (see below) to remove primer-dimers or small DNA molecules (<300 bp).
- 30% PEG 8000/30 mM MgCl<sub>2</sub> Solution is provided to purify PCR products away from other DNA <300 bp in size, including primer-dimers. Run a 25 µl PCR reaction, dilute the PCR reaction 4-fold with TE [10 mM Tris-HCl (pH 7.5-8), 1 mM EDTA], add 1/2 volume of 30% PEG 8000/30 mM MgCl<sub>2</sub> Solution (final concentrations of 10% PEG, 10 mM MgCl<sub>2</sub>), and vortex. Centrifuge 15 minutes at full speed in a microcentrifuge. Carefully remove supernatant and suspend the clear pellet in TE to >10 ng/µl.

**Important:** Use the recommended proportion of PEG/MgCl<sub>2</sub> to ensure that correct-sized products are removed.

- For BP reactions, the most efficient substrates are **linear** *attB* products (PCR products or expression clones) and supercoiled *attP*-containing donor vectors. Supercoiled or relaxed *attB* substrates may be used but will react less efficiently than linear *attB* substrates.
- To increase the number of colonies containing the desired entry clone, increase the incubation time from the recommended 1 hour to 4-6 hours (typically 2-3 fold more colonies) or overnight (typically 5-10 fold more colonies). Longer incubations are recommended for genes ≥ 5 kb to increase the yield of colonies.
- We recommend using 20-50 fmol of PCR product per 10 µl reaction (where a 1 kb PCR product is ~0.65 ng/fmol). Increasing the amount of PCR product generally yields more colonies; however, do not exceed ~250 ng of PCR product per 10 µl reaction.

## Procedures

### BP Reaction

BP Clonase™ II enzyme mix is supplied as a 5X solution. If you wish to scale the reaction volume, make sure the BP Clonase™ II enzyme mix is at a final concentration of 1X. For a positive control, use 100 ng (2 µl) of pEXP7-tet.

1. Add the following components to a 1.5 ml microcentrifuge tube at room temperature and mix:

<i>attB</i> -PCR product (≥10 ng/µl; final amount ~15-150 ng)	1-7 µl
Donor vector (150 ng/µl)	1 µl
TE buffer, pH 8.0	to 8 µl
2. Thaw on ice the BP Clonase™ II enzyme mix for about 2 minutes. Vortex the BP Clonase™ II enzyme mix briefly twice (2 seconds each time).
3. To each sample (Step 1, above), add 2 µl of BP Clonase™ II enzyme mix to the reaction and mix well by vortexing briefly twice. Microcentrifuge briefly.
4. Return BP Clonase™ II enzyme mix to -20°C or -80°C storage.
5. Incubate reactions at 25°C for 1 hour.
6. Add 1 µl of the Proteinase K solution to each sample to terminate the reaction. Vortex briefly. Incubate samples at 37°C for 10 minutes.

### Transformation

1. Transform 1 µl of each BP reaction into 50 µl of One Shot® OmniMAX™ 2 T1 Phage-Resistant Cells (Catalog no. C8540-03). Incubate on ice for 30 minutes. Heat-shock cells by incubating at 42°C for 30 seconds. Add 250 µl of S.O.C. Medium and incubate at 37°C for 1 hour with shaking. Plate 20 µl and 100 µl of each transformation onto selective plates. **Note:** Any competent cells with a transformation efficiency of  $>1.0 \times 10^8$  transformants/µg may be used.
2. Transform 1 µl of pUC19 DNA (10 ng/ml) into 50 µl of One Shot® OmniMAX™ 2 T1 Phage-Resistant Cells as described above. Plate 20 µl and 100 µl on LB plates containing 100 µg/ml ampicillin.

### Expected Results

An efficient BP recombination reaction will produce >1500 colonies if the entire BP reaction is transformed and plated.

## References

1. Landy, A. (1989) *Ann. Rev. Biochem.* 58, 913.
2. Bernard, P. and Couturier, M. (1992) *J. Mol. Biol.* 226, 735.
3. Miki, T., Park, J.A., Nagao, K., Murayama, N., and Horiuchi, T. (1992) *J. Mol. Biol.* 225, 39.

## Limited Use Label License No. 19: Gateway<sup>®</sup> Cloning Products

This product and its use is the subject of one or more of U.S. Patent Nos. 5,888,732, 6,143,557, 6,171,861, 6,270,969, 6,277,608, and 6,720,140 and/or other pending U.S. and foreign patent applications owned by Invitrogen Corporation. The purchase of this product conveys to the buyer the non-transferable right to use the purchased amount of the product and components of the product in research conducted by the buyer (whether the buyer is an academic or for profit entity). The purchase of this product does not convey a license under any method claims in the foregoing patents or patent applications, or to use this product with any recombination sites other than those purchased from Invitrogen Corporation or its authorized distributor. The right to use methods claimed in the foregoing patents or patent applications with this product for research purposes only can only be acquired by the use of Clonase<sup>™</sup> purchased from Invitrogen Corporation or its authorized distributors. The buyer cannot modify the recombination sequence(s) contained in this product for any purpose. The buyer cannot sell or otherwise transfer (a) this product, (b) its components, or (c) materials made by the employment of this product or its components to a third party or otherwise use this product or its components or materials made by the employment of this product or its components for Commercial Purposes. The buyer may transfer information or materials made through the employment of this product to a scientific collaborator, provided that such transfer is not for any Commercial Purpose, and that such collaborator agrees in writing (a) not to transfer such materials to any third party, and (b) to use such transferred materials and/or information solely for research and not for Commercial Purposes. Notwithstanding the preceding, any buyer who is employed in an academic or government institution may transfer materials made with this product to a third party who has a license from Invitrogen under the patents identified above to distribute such materials. Transfer of such materials and/or information to collaborators does not convey rights to practice any methods claimed in the foregoing patents or patent applications. Commercial Purposes means any activity by a party for consideration and may include, but is not limited to: (1) use of the product or its components in manufacturing; (2) use of the product or its components to provide a service, information, or data; (3) use of the product or its components for therapeutic, diagnostic or prophylactic purposes; or (4) resale of the product or its components, whether or not such product or its components are resold for use in research. Invitrogen Corporation will not assert a claim against the buyer of infringement of the above patents based upon the manufacture, use or sale of a therapeutic, clinical diagnostic, vaccine or prophylactic product developed in research by the buyer in which this product or its components was employed, provided that none of (i) this product, (ii) any of its components, or (iii) a method claim of the foregoing patents, was used in the manufacture of such product. Invitrogen Corporation will not assert a claim against the buyer of infringement of the above patents based upon the use of this product to manufacture a protein for sale, provided that no method claim in the above patents was used in the manufacture of such protein. If the purchaser is not willing to accept the limitations of this limited use statement, Invitrogen is willing to accept return of the product with a full refund. For information on purchasing a license to use this product for purposes other than those permitted above, contact Licensing Department, Invitrogen Corporation, 1600 Faraday Avenue, Carlsbad, California 92008. Phone (760) 603-7200.

## Limited Use Label License No. 28: CMV Promoter

The use of the CMV promoter is covered under U.S. Patent Nos. 5,168,062 and 5,385,839 owned and licensed by the University of Iowa Research Foundation and is sold for research use only. Commercial users must obtain a license to these patents directly from the University of Iowa Research Foundation (UIRF), 214 Technology Innovation Center, Iowa City, Iowa 52242. For further information, please contact the Associate Director of UIRF, at 319-335-4546.

©2004-2005 Invitrogen Corporation. All rights reserved.