

June 2004

# MinElute<sup>®</sup> Handbook

For

MinElute PCR Purification Kit

MinElute Gel Extraction Kit

MinElute Reaction Cleanup Kit



WWW.QIAGEN.COM

Trademarks: QIAGEN®, MinElute® (QIAGEN Group), Tween® (ICI Americas Inc.). The PCR process is covered by U.S. patents 4,683,195 and 4,683,202 and foreign equivalents owned by Hoffmann-La Roche AG.

© 2004 QIAGEN, all rights reserved.

# Contents

<b>Kit Contents</b>	<b>4</b>
<b>Storage Conditions</b>	<b>4</b>
<b>Product Use Limitations</b>	<b>5</b>
<b>Product Warranty and Satisfaction Guarantee</b>	<b>5</b>
<b>Quality Control</b>	<b>5</b>
<b>Technical Assistance</b>	<b>5</b>
<b>Safety Information</b>	<b>6</b>
<b>Introduction</b>	<b>7</b>
<b>Specifications</b>	<b>8</b>
<b>The MinElute Principle</b>	<b>10</b>
<b>Protocols</b>	<b>16</b>
<b>MinElute PCR Purification Kit Protocols</b>	<b>16</b>
■ using a microcentrifuge	16
■ using a vacuum manifold	17
<b>MinElute Gel Extraction Kit Protocols</b>	<b>19</b>
■ using a microcentrifuge	19
■ using a vacuum manifold	21
<b>MinElute Reaction Cleanup Kit Protocols</b>	<b>24</b>
■ using a microcentrifuge	24
■ using a vacuum manifold	26
<b>Troubleshooting Guide</b>	<b>28</b>
<b>Appendix: QIAvac Vacuum Manifolds</b>	<b>30</b>
<b>Handling Guidelines for QIAvac 24</b>	<b>30</b>
<b>Handling Guidelines for QIAvac 6S</b>	<b>31</b>
<b>References</b>	<b>32</b>
<b>Ordering Information</b>	<b>33</b>
<b>QIAGEN Companies and Distributors</b>	<b>35</b>

## Kit Contents

<b>MinElute® PCR Purification Kits</b>	<b>(50)</b>	<b>(250)</b>
<b>Catalog no.</b>	<b>28004</b>	<b>28006</b>
MinElute Spin Columns	50	250
Buffer PB*	30 ml	150 ml
Buffer PE (concentrate)	2 x 6 ml	55 ml
Buffer EB	15 ml	15 ml
Collection Tubes (2 ml)	50	250
Handbook	1	1
<b>MinElute Gel Extraction Kits</b>	<b>(50)</b>	<b>(250)</b>
<b>Catalog no.</b>	<b>28604</b>	<b>28606</b>
MinElute Spin Columns	50	250
Buffer QG*	2 x 50 ml	2 x 250 ml
Buffer PE (concentrate)	2 x 6 ml	55 ml
Buffer EB	15 ml	15 ml
Collection Tubes (2 ml)	50	250
Handbook	1	1
<b>MinElute Reaction Cleanup Kits</b>	<b>(50)</b>	<b>(250)</b>
<b>Catalog no.</b>	<b>28204</b>	<b>28206</b>
MinElute Spin Columns	50	250
Buffer ERC*	15 ml	85 ml
Buffer PE	2 x 6 ml	55 ml
Buffer EB	15 ml	15 ml
Collection Tubes (2 ml)	50	250
Handbook	1	1

\* Buffers PB, QG, and ERC contain chaotropic salts which are irritants. Take appropriate laboratory safety measures and wear gloves when handling.

## Storage Conditions

Upon arrival, open the kit and store MinElute spin columns at 2–8°C. The remaining kit components can be stored at room temperature (15–25°C). Under these conditions, MinElute Kits can be stored for up to 12 months without showing any reduction in performance and quality. The entire kit can be stored at 2–8°C, but in this case buffers should be redissolved before use. Make sure that all buffers are at room temperature when used.

## Product Use Limitations

MinElute Kits are developed, designed and sold for research purposes only. They are not to be used for human diagnostic or drug purposes or to be administered to humans unless expressly cleared for that purpose by the Food and Drug Administration in the USA or the appropriate regulatory authorities in the country of use. All due care and attention should be exercised in the handling of many of the materials described in this text.

## Product Warranty and Satisfaction Guarantee

QIAGEN guarantees the performance of all products in the manner described in our product literature. The purchaser must determine the suitability of the product for its particular use. Should any product fail to perform satisfactorily due to any reason other than misuse, QIAGEN will replace it free of charge or refund the purchase price. We reserve the right to change, alter, or modify any product to enhance its performance and design. If a QIAGEN product does not meet your expectations, simply call your local Technical Service Department. We will credit your account or exchange the product — as you wish.

A copy of QIAGEN terms and conditions can be obtained on request, and is also provided on the back of our invoices. If you have questions about product specifications or performance, please call QIAGEN Technical Services or your local distributor.

## Quality Control

As part of the stringent QIAGEN quality assurance program, the performance of MinElute Kits is monitored routinely and on a lot-to-lot basis. MinElute Kits are tested by isolation of DNA fragments of various sizes from either aqueous solution or agarose gel. The quality of the isolated DNA is checked by several assays commonly used for nucleic acids. The DNA binding capacity of MinElute spin columns is tested by determining the recovery from a specific amount of loaded DNA.

## Technical Assistance

At QIAGEN we pride ourselves on the quality and availability of our technical support. Our Technical Service Departments are staffed by experienced scientists with extensive practical and theoretical expertise in molecular biology and the use of QIAGEN products. If you have any questions or experience any problems regarding any aspect of MinElute Kits, or QIAGEN products in general, please do not hesitate to contact us.

QIAGEN customers are a major source of information regarding advanced or specialized uses of our products. This information is helpful to other scientists as well as to the researchers at QIAGEN. We therefore encourage you to contact us if you have any suggestions about product performance or new applications and techniques.

For technical assistance and more information please call one of the QIAGEN Technical Service Departments or local distributors (see back cover).

## Safety Information

When working with chemicals, always wear a suitable lab coat, disposable gloves, and protective goggles. For more information, please consult the appropriate material safety data sheets (MSDSs). These are available online in convenient and compact PDF format at [www.qiagen.com/ts/msds.asp](http://www.qiagen.com/ts/msds.asp) where you can find, view, and print the MSDS for each QIAGEN kit and kit component.

**CAUTION: DO NOT add bleach or acidic solutions directly to the sample-preparation waste.**

Buffers PB and ERC contain guanidine hydrochloride, which can form highly reactive compounds when combined with bleach.

In case liquid containing these buffers is spilled, clean with suitable laboratory detergent and water. If the spilled liquid contains potentially infectious agents, clean the affected area first with laboratory detergent and water, and then with 1% (v/v) sodium hypochlorite.

The following risk and safety phrases apply to the components of the MinElute DNA Cleanup System

### Buffers PB

Contains guanidine hydrochloride and isopropanol: harmful, irritant, flammable. Risk and safety phrases: \* R10-22-36/38, S13-23-26-36/37/39-46

### Buffer ERC

Contains guanidine hydrochloride and isopropanol: harmful, irritant, flammable. Risk and safety phrases: \* R10-22-36/38, S23-26-36/37/39-46

### Buffer QG

Contains guanidine thiocyanate: harmful, irritant. Risk and safety phrases: \* R20/21/22-32, S13-26-36-46

### 24-hour emergency information

Emergency medical information in English, French, and German can be obtained 24 hours a day from:

Poison Information Center Mainz, Germany

Tel: +49-6131-19240

\* R10: Flammable. R22: Harmful if swallowed. R20/21/22: Harmful by inhalation, in contact with skin and if swallowed. R32: Contact with acids liberates very toxic gas. R36/38: Irritating to eyes and skin. S13: Keep away from food, drink and animal feedingstuffs. S23: Do not breathe vapour/spray. S26: In case of contact with eyes, rinse immediately with plenty of water and seek medical advice. S36: Wear suitable protective clothing. S36/37/39: Wear suitable protective clothing, gloves and eye/face protection. S46: If swallowed, seek medical advice immediately and show the container or label.

## Introduction

The MinElute DNA Cleanup system is specially designed for fast and easy isolation of DNA fragments from PCR reactions, agarose gels, or enzymatic reactions with an extremely small elution volume of only 10  $\mu$ l. Designed for rapid DNA cleanup, with extremely high end-concentration of DNA, the MinElute system includes:

- **MinElute PCR Purification Kits** for direct purification of double-stranded PCR products (70 bp – 4 kb) from amplification reactions.
- **MinElute Gel Extraction Kits** for extraction of DNA fragments (70 bp – 4 kb) from standard or low-melt agarose gels in TAE (Tris-acetate/EDTA) or TBE (Tris-borate/EDTA) buffer.
- **MinElute Reaction Cleanup Kits** to purify double-stranded DNA fragments (70 bp – 4 kb) from all enzymatic reactions.

The MinElute Reaction Cleanup Kit can be used for DNA cleanup from all enzymatic reactions, including:

- Dephosphorylation
- Restriction enzyme digestion
- Ligation
- Primed synthesis
- Endlabeling
- Nick translation

During DNA cleanup using MinElute Reaction Cleanup Kits, all enzymes are removed, independent of size and secondary structure. Table 1 lists examples of enzymes that are used in the above reactions and are removed, together with salts and oligomers, during the MinElute Reaction Cleanup procedure.

**Table 1. Enzymes Commonly Used in Enzymatic Manipulation of DNA**

Protein	Molecular weight per enzyme subunit (kDa)
DNA Polymerase I	109
Klenow fragment	62
Calf intestinal alkaline phosphatase (CIP)	69
T4 DNA ligase	55
T4 Polynucleotide kinase	35
Terminal transferase	32
DNase I	31

## Specifications

	MinElute PCR Purification Kit	MinElute Gel Extraction Kit	MinElute Reaction Cleanup Kit
Maximum binding capacity:	5 µg	5 µg	5 µg
Recovery of DNA:	80% (70 bp – 4kb)	80% (70 bp – 4 kb)	80% (70 bp – 4 kb)
Maximum weight of gel slice:	—	400 mg	—
Elution volume:	10 µl	10 µl	10 µl
Volume of eluate:	9 µl	9 µl	9 µl
Capacity of column reservoir:	800 µl	800 µl	800 µl
<b>Recovered:</b>			
dsDNA	70 bp – 4 kb	70 bp – 4 kb	70 bp – 4 kb
<b>Removed:</b>			
<40mers	<b>YES</b>	n.a.	<b>YES</b>


n.a.: not applicable

MinElute Kits provide high yields of pure nucleic acids in minimal elution volumes. The highly concentrated DNA is suitable for direct use in applications such as:

- Ligation and transformation
- Radioactive and fluorescent sequencing
- Restriction enzyme digestion
- Labeling
- In vitro transcription
- Amplification
- Microinjection
- Hybridization



## DNA Fragment Binding-Size Range

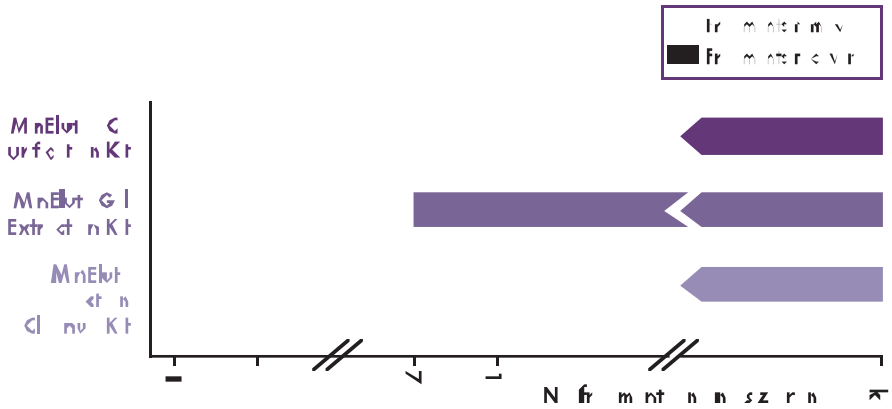


Recoveries of DNA fragments in the size range between “removed” and “recovered” are not defined.

## The MinElute Principle

The MinElute system combines the convenience of spin-column technology with the selective binding properties of a uniquely designed silica-gel membrane. MinElute columns are designed to give high end-concentrations of purified DNA fragments for subsequent reactions. Special buffers provided with each kit are optimized for efficient recovery of DNA and removal of contaminants in each specific application. DNA adsorbs to the silica-membrane in the presence of high salt while contaminants pass through the column. Impurities are efficiently washed away, and the pure DNA is eluted with Tris buffer\* or water (see page 15). The MinElute columns offer two handling options — as an alternative to processing the spin columns in a microcentrifuge, they can also be used on any commercial vacuum manifold with luer connectors, e.g., QIAvac 24 (or QIAvac 6S with QIAvac Luer Adapters). The uniquely designed MinElute column with specialized membrane assembly allows purification of concentrated DNA fragments in high yields, in as little as 9  $\mu$ l eluate (Figure 1).

### Unique Membrane Assembly Enables Low Elution Volumes

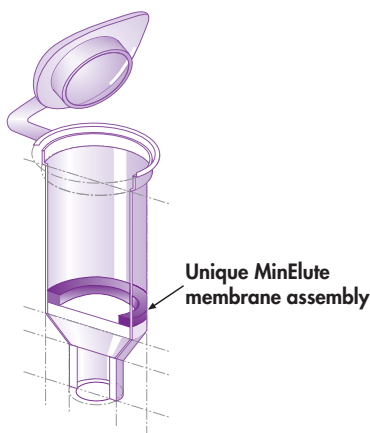


**Figure 1** Cross-section of MinElute column (utility model pending).

\* When working with chemicals, always wear a suitable lab coat, disposable gloves, and protective goggles. For more information, please consult the appropriate material safety data sheets (MSDSs) available from the product supplier.

## Adsorption to MinElute membrane — salt and pH dependence

The MinElute silica-gel membrane is uniquely adapted to isolate DNA from both aqueous solutions and agarose gels, and up to 5 µg DNA can bind to each MinElute column. The binding buffers in MinElute Spin Kits provide the correct salt concentration and pH for adsorption of DNA to the MinElute membrane. The adsorption of nucleic acids to silica-gel surfaces occurs only in the presence of a high concentration of chaotropic salts (1), which modify the structure of water (2). Adsorption of DNA to silica also depends on pH. Adsorption is typically 80% if the pH is  $\leq 7.5$ , and is reduced drastically at higher pH (Figure 2). If the loading mixture pH is  $>7.5$ , the optimal pH for DNA binding can be obtained by adding a small volume of 3 M sodium acetate\*, pH 5.0.

### pH Dependence of DNA Adsorption to MinElute Membranes

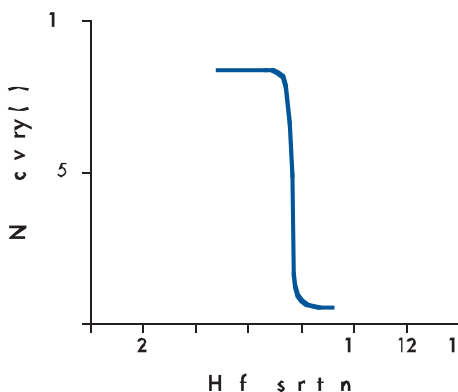

**Figure 2** pH dependence of DNA adsorption to silica. 1 µg of a 2.9 kb DNA fragment was adsorbed at different pHs and eluted with Buffer EB (10 mM Tris-Cl, pH 8.5). The graph shows the percentage of DNA recovery, reflecting the relative adsorption efficiency, versus pH of adsorption.

### Optimized binding buffers for every DNA cleanup task

MinElute Kits contain identical MinElute columns but different binding buffers optimized for each specific application:

- Buffer PB in the MinElute PCR Purification Kit allows the efficient binding of double-stranded PCR products as small as 70 bp and the removal of primers up to 40 nucleotides.
- Buffer QG in the MinElute Gel Extraction Kit solubilizes the agarose gel slice and provides the appropriate conditions for binding of DNA to the silica membrane.
- Buffer ERC in the MinElute Reaction Cleanup Kit allows the efficient binding of double stranded DNA as small as 70 bp and the removal of enzymes, salts, and oligomers.
- Buffer PB and Buffer QG are also available separately (see Ordering Information).

## pH indicator in Buffers QG and ERC

The binding and solubilization buffer QG and binding buffer ERC are specially optimized for use with the MinElute silica-gel membrane. The buffers contain a pH indicator, allowing easy determination of the optimal pH for DNA binding. DNA adsorption requires a pH  $\leq 7.5$ , and the pH indicator in Buffers QG and ERC appears yellow in this range. If the pH is  $>7.5$ , which can occur if during agarose gel electrophoresis, the electrophoresis buffer had been used repeatedly or incorrectly prepared, or if the buffer used in an enzymatic reaction is strongly basic and has a high buffering capacity, the binding mixture turns orange or violet (Figure 3). This means that the pH of the sample exceeds the buffering capacity of Buffer QG or ERC and DNA adsorption will be inefficient. In these cases, the pH of the binding mixture can easily be corrected by addition of a small volume of 3 M sodium acetate\*, pH 5.0, before proceeding with the protocol. In addition, in the MinElute Gel Extraction Kit procedure, the color of the binding mixture allows easy visualization of any unsolubilized agarose, ensuring complete solubilization and maximum yields. The indicator dye does not interfere with DNA binding and is completely removed during the cleanup procedure. Buffers QG and ERC do not contain sodium iodide (NaI). Residual NaI may be difficult to remove from DNA samples, and reduces the efficiency of subsequent enzymatic reactions such as blunt-end ligations.

### Indicator Enables Easy Checking of the Optimal pH



**Figure 3** Indicator dye in solubilization and binding Buffers QG and ERC identifies optimal pH for DNA binding.

\* When working with chemicals, always wear a suitable lab coat, disposable gloves, and protective goggles. For more information, please consult the appropriate material safety data sheets (MSDSs) available from the product supplier.

## Washing

During the DNA adsorption step, unwanted primers and impurities, such as salts, enzymes, unincorporated nucleotides, agarose, dyes, ethidium bromide, DMSO, oils, and detergents (e.g., Tween® 20) do not bind to the silica membrane, but flow through the column. Salts are quantitatively washed away by the ethanol-containing Buffer PE. Any residual Buffer PE, which may interfere with subsequent enzymatic reactions, is removed by an additional centrifugation step.

## Elution in low-salt solutions

Elution efficiency is strongly dependent on the salt concentration and pH of the elution buffer. Contrary to adsorption, elution is most efficient under basic conditions and low salt concentrations. DNA is eluted with 10 µl of the provided Buffer EB (10 mM Tris-Cl, pH 8.5), or water. The volume of the final eluate is 9 µl. The maximum elution efficiency is achieved between pH 7.0 and 8.5. When using water to elute, make sure that the pH is within this range. In addition, DNA must be stored at -20°C when eluted with water since DNA may degrade in the absence of a buffering agent. Elution with TE (10 mM Tris-Cl, 1 mM EDTA, pH 8.0) is possible, but not recommended because EDTA may inhibit subsequent enzymatic reactions.

## Agarose gel analysis of yield

Yields of DNA following cleanup can be analyzed by agarose gel analysis. Table 2 shows the total yield obtained following extraction of 1 µg or 0.5 µg starting DNA from an agarose gel with a recovery of 80% using the MinElute Gel Extraction Kit. The corresponding amount of DNA in a 1 µl aliquot from 10 µl elution volume (9 µl eluate) is indicated.

**Table 2. Amount of DNA in 1 µl Aliquots of a 9 µl Eluate Following MinElute Purification**

Starting DNA	Recovery	Total yield (10 µl elution volume)	Amount of DNA in 1 µl (9 µl eluate)
1 µg	80%	800 ng	89 ng
0.5 µg	80%	400 ng	44 ng

## Quantitation of DNA fragments

DNA fragments can be quantitated by running a sample alongside standards containing known quantities of the same-sized DNA fragment. The amount of sample DNA loaded can be estimated by visual comparison of the band intensity with that of the standards (Figure 4).




**Figure 4** An unknown amount of a 5.5 kb DNA fragment (**U**) was run alongside known quantities (as indicated in ng) of the same DNA fragment. The unknown sample contained 75–100 ng DNA, as estimated by visual comparison with the standards. **M**: 1 kb DNA ladder.

## Applications using MinElute purified DNA

DNA purified by the MinElute System is significantly more concentrated than DNA purified by other methods. The highly concentrated DNA makes small reaction volumes possible for downstream applications, leading to increased efficiency (e.g., in ligations). DNA purified with MinElute Kits is suitable for any subsequent application, such as ligation and transformation, radioactive and fluorescent sequencing, restriction enzyme digestion, labeling, hybridization, PCR, in vitro transcription, or microinjection.

## The MinElute Purification Procedure

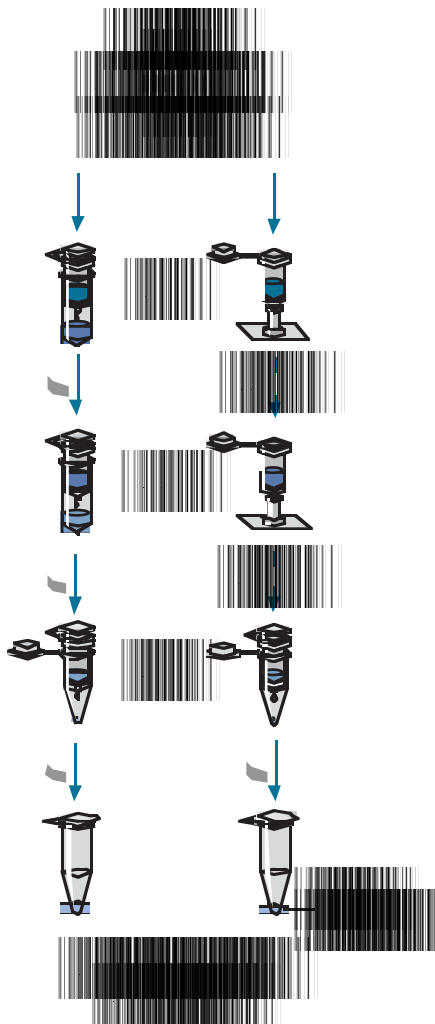

# MinElute PCR Purification Kit Protocol

## using a microcentrifuge

This protocol is designed to purify double-stranded DNA fragments from PCR reactions resulting in high end-concentrations of DNA (see page 10). Fragments ranging from 70 bp to 4 kb are purified from primers, nucleotides, polymerases, and salts using MinElute spin columns in a microcentrifuge.

### Important points before starting

- Add ethanol\* (96–100%) to Buffer PE before use (see bottle label for volume).
- All centrifuge steps are at  $\geq 10,000 \times g$  (~13,000 rpm) in a conventional tabletop microcentrifuge.

1. **Add 5 volumes of Buffer PB to 1 volume of the PCR reaction and mix. It is not necessary to remove mineral oil or kerosene.**

For example, add 250  $\mu$ l of Buffer PB to 50  $\mu$ l PCR reaction (not including oil).

2. **Place a MinElute column in a provided 2 ml collection tube in a suitable rack.**
3. **To bind DNA, apply the sample to the MinElute column and centrifuge for 1 min.**  
For maximum recovery, transfer all traces of sample to the column.
4. **Discard flow-through. Place the MinElute column back into the same tube.**
5. **To wash, add 750  $\mu$ l Buffer PE to the MinElute column and centrifuge for 1 min.**
6. **Discard flow-through and place the MinElute column back in the same tube. Centrifuge the column for an additional 1 min at maximum speed.**

**IMPORTANT:** Residual ethanol from Buffer PE will not be completely removed unless the flow-through is discarded before this additional centrifugation.

7. **Place the MinElute column in a clean 1.5 ml microcentrifuge tube.**
8. **To elute DNA, add 10  $\mu$ l Buffer EB (10 mM Tris-Cl, pH 8.5) or H<sub>2</sub>O to the center of the membrane, let the column stand for 1 min, and then centrifuge for 1 min.**

**IMPORTANT:** Ensure that the elution buffer is dispensed directly onto the center of the membrane for complete elution of bound DNA. The average eluate volume is 9  $\mu$ l from 10  $\mu$ l elution buffer volume.

Elution efficiency is dependent on pH. The maximum elution efficiency is achieved between pH 7.0 and 8.5. When using water, make sure that the pH value is within this range, and store DNA at  $-20^{\circ}\text{C}$  as DNA may degrade in the absence of a buffering agent. The purified DNA can also be eluted in TE\* (10 mM Tris-Cl, 1 mM EDTA, pH 8.0), but the EDTA may inhibit subsequent enzymatic reactions.

\* When working with chemicals, always wear a suitable lab coat, disposable gloves, and protective goggles. For more information, please consult the appropriate material safety data sheets (MSDSs) available from the product supplier.



# MinElute PCR Purification Kit Protocol

## using a vacuum manifold

MinElute spin columns can also be used on any vacuum manifold with luer connectors, e.g., QIAvac 24 (or QIAvac 6S with Luer Adapters). The following protocol is designed to purify double-stranded DNA fragments from PCR reactions resulting in high end-concentrations of DNA (see page 10). Fragments ranging from 70 bp to 4 kb are purified from primers, nucleotides, polymerases and salts using vacuum-driven sample processing.

### Important points before starting

- Add ethanol (96–100%) to Buffer PE before use (see bottle label for volume).
- Switch off vacuum between steps to ensure that a consistent, even vacuum is applied during manipulations.

### 1. Add 5 volumes of Buffer PB to 1 volume of the PCR reaction and mix. It is not necessary to remove mineral oil or kerosene.

For example, add 250  $\mu$ l of Buffer PB to 50  $\mu$ l PCR reaction (not including oil).

### 2. Prepare the vacuum manifold and MinElute columns:

#### 2a. QIAvac 24 (see page 30):

Place the QIAvac 24 lid on top of the QIAvac 24 base. Make sure that the gasket fits tightly in the groove of the QIAvac 24 lid.

Insert up to 24 MinElute spin columns into the luer extensions of the QIAvac 24. Close unused positions with luer caps and connect QIAvac 24 to a vacuum source.

#### 2b. QIAvac 6S manifold (see page 31):

Open QIAvac 6S lid. Place QIAvac Luer Adapter(s), or blanks to seal unused slots, into the slots of QIAvac top plate, and close the QIAvac 6S lid. Place the waste tray inside the QIAvac base, and place the top plate squarely over the base. Attach the QIAvac 6S to a vacuum source.

Insert each MinElute column into a luer connector on the Luer Adapter(s) in the manifold. Seal unused luer connectors with plugs provided with the QIAvac Luer Adapter Set.

#### 2c. Other vacuum manifolds: follow the supplier's instructions. Insert each MinElute column into a luer connector.

### 3. To bind DNA, load the samples into the MinElute columns by decanting or pipetting, and apply vacuum. After the samples have passed through the columns, switch off the vacuum source.

For maximum recovery, transfer all traces of sample to the column.

The maximum loading volume of the column is 800  $\mu$ l. For sample volumes greater than 800  $\mu$ l simply load again.

4. To wash, add 750  $\mu$ l of Buffer PE to each MinElute column and apply vacuum.
5. Transfer each MinElute column to a microcentrifuge tube or the provided 2 ml collection tubes. Centrifuge for 1 min at  $\geq 10,000 \times g$  (~13,000 rpm).

**IMPORTANT:** This spin is necessary to remove residual ethanol (Buffer PE).

6. Place each MinElute column into a clean 1.5 ml microcentrifuge tube.
7. To elute DNA, add 10  $\mu$ l of Buffer EB (10 mM Tris-Cl, pH 8.5) or H<sub>2</sub>O to the center of each membrane, let the columns stand for 1 min, and then centrifuge.

**IMPORTANT:** Ensure that the elution buffer is dispensed directly onto the center of the membrane for complete elution of bound DNA. The average eluate volume is 9  $\mu$ l from 10  $\mu$ l elution buffer volume.

Elution efficiency is dependent on pH. The maximum elution efficiency is achieved between pH 7.0 and 8.5. When using water, make sure that the pH value is within this range, and store DNA at  $-20^{\circ}\text{C}$  as DNA may degrade in the absence of a buffering agent. The purified DNA can also be eluted in TE\* (10 mM Tris-Cl, 1 mM EDTA, pH 8.0), but the EDTA may inhibit subsequent enzymatic reactions.

\* When working with chemicals, always wear a suitable lab coat, disposable gloves, and protective goggles. For more information, please consult the appropriate material safety data sheets (MSDSs) available from the product supplier.

# MinElute Gel Extraction Kit Protocol

## using a microcentrifuge

This protocol is designed to extract and purify DNA of 70 bp to 4 kb from standard or low-melt agarose gels in TAE or TBE buffer resulting in high end-concentrations of DNA. Up to 400 mg agarose can be processed per MinElute column.

### Important points before starting

- The yellow color of Buffer QG indicates a pH  $\leq 7.5$ .
- Add ethanol\* (96–100%) to Buffer PE before use (see bottle label for volume).
- Isopropanol\* (100%) and a heating block or water bath at 50°C are required.
- All centrifugation steps are carried out at  $\geq 10,000 \times g$  (~13,000 rpm) in a conventional table-top microcentrifuge.
- 3 M sodium acetate\*, pH 5.0, may be required.

### 1. Excise the DNA fragment from the agarose gel with a clean, sharp scalpel.

Minimize the size of the gel slice by removing extra agarose.

### 2. Weigh the gel slice in a colorless tube. Add 3 volumes of Buffer QG to 1 volume of gel (100 mg ~ 100 $\mu$ l).

For example, add 300  $\mu$ l of Buffer QG to each 100 mg of gel. For >2% agarose gels, add 6 volumes of Buffer QG. The maximum amount of gel slice per spin column is 400 mg; for gel slices >400 mg use more than one MinElute column.

### 3. Incubate at 50°C for 10 min (or until the gel slice has completely dissolved). To help dissolve gel, mix by vortexing the tube every 2–3 min during the incubation.

**IMPORTANT:** Solubilize agarose completely. For >2% gels, increase incubation time.

### 4. After the gel slice has dissolved completely, check that the color of the mixture is yellow (similar to Buffer QG without dissolved agarose).

**Note:** If the color of the mixture is orange or violet, add 10  $\mu$ l of 3 M sodium acetate, pH 5.0, and mix. The color of the mixture will turn to yellow.

The adsorption of DNA to the membrane is efficient only at pH  $\leq 7.5$ . Buffer QG contains a pH indicator which is yellow at pH  $\leq 7.5$  and orange or violet at higher pH, allowing easy determination of the optimal pH for DNA binding.

### 5. Add 1 gel volume of isopropanol to the sample and mix by inverting the tube several times.

For example, if the agarose gel slice is 100 mg, add 100  $\mu$ l isopropanol. Do not centrifuge the sample at this stage.

\* When working with chemicals, always wear a suitable lab coat, disposable gloves, and protective goggles. For more information, please consult the appropriate material safety data sheets (MSDSs) available from the product supplier.

6. Place a MinElute column in a provided 2 ml collection tube in a suitable rack.
7. To bind DNA, apply the sample to the MinElute column, and centrifuge for 1 min.

For maximum recovery, transfer all traces of sample to the column. The maximum volume of the column reservoir is 800  $\mu$ l. For sample volumes of more than 800  $\mu$ l, simply load and spin again.

8. Discard the flow-through and place the MinElute column back in the same collection tube.
9. Add 500  $\mu$ l of Buffer QG to the spin column and centrifuge for 1 min.
10. Discard the flow-through and place the MinElute column back in the same collection tube.
11. To wash, add 750  $\mu$ l of Buffer PE to the MinElute column and centrifuge for 1 min.

**Note:** If the DNA will be used for salt sensitive applications, such as blunt-end ligation and direct sequencing, let the column stand 2–5 min after addition of Buffer PE, before centrifuging.

12. Discard the flow-through and centrifuge the MinElute column for an additional 1 min at  $\geq 10,000 \times g$  (~13,000 rpm).

**IMPORTANT:** Residual ethanol from Buffer PE will not be completely removed unless the flow-through is discarded before this additional centrifugation.

13. Place the MinElute column into a clean 1.5 ml microcentrifuge tube.
14. To elute DNA, add 10  $\mu$ l of Buffer EB (10 mM Tris-Cl, pH 8.5) or H<sub>2</sub>O to the center of the membrane, let the column stand for 1 min, and then centrifuge for 1 min.

**IMPORTANT:** Ensure that the elution buffer is dispensed directly onto the center of the membrane for complete elution of bound DNA. The average eluate volume is 9  $\mu$ l from 10  $\mu$ l elution buffer volume.

Elution efficiency is dependent on pH. The maximum elution efficiency is achieved between pH 7.0 and 8.5. When using water, make sure that the pH value is within this range, and store DNA at  $-20^{\circ}\text{C}$  as DNA may degrade in the absence of a buffering agent. The purified DNA can also be eluted in TE\* (10 mM Tris-Cl, 1 mM EDTA, pH 8.0), but the EDTA may inhibit subsequent enzymatic reactions.

\* When working with chemicals, always wear a suitable lab coat, disposable gloves, and protective goggles. For more information, please consult the appropriate material safety data sheets (MSDSs) available from the product supplier.

# MinElute Gel Extraction Kit Protocol

## using a vacuum manifold

MinElute columns can also be used on any vacuum manifold with luer connectors, e.g., QIAvac 24 (or QIAvac 6S with Luer Adapters). The following protocol is designed to extract and purify DNA of 70 bp to 4 kb from standard or low-melt agarose gels in TAE or TBE buffer using vacuum-driven processing resulting in high end-concentrations of DNA. Up to 400 mg agarose can be processed per MinElute column.

### Important points before starting

- The yellow color of Buffer QG indicates a pH  $\leq 7.5$ .
- Add ethanol\* (96–100%) to Buffer PE before use (see bottle label for volume).
- Isopropanol\* (100%) and a heating block or water bath at 50°C are required.
- Switch off vacuum between steps to ensure that a consistent, even vacuum is applied during manipulations.
- 3 M sodium acetate\*, pH 5.0, may be required.

### 1. Excise the DNA fragment from the agarose gel with a clean, sharp scalpel.

Minimize the size of the gel slice by removing extra agarose.

### 2. Weigh the gel slice in a colorless tube. Add 3 volumes of Buffer QG to 1 volume of gel (100 mg ~ 100 $\mu$ l).

For example, add 300  $\mu$ l of Buffer QG to each 100 mg of gel. For >2% agarose gels, add 6 volumes of Buffer QG. The maximum amount of gel slice per spin column is 400 mg; for gel slices >400 mg use more than one spin column.

### 3. Incubate at 50°C for 10 min (or until the gel slice has completely dissolved). To help dissolve gel, mix by vortexing the tube every 2–3 min during the incubation.

**IMPORTANT:** Solubilize agarose completely. For >2% gels, increase incubation time.

### 3a. During the incubation, prepare the vacuum manifold and MinElute columns:

#### QIAvac 24 (see page 30):

Place the QIAvac 24 lid on top of the QIAvac 24 base. Make sure that the gasket fits tightly in the groove of the QIAvac 24 lid.

Insert up to 24 MinElute columns into the luer extensions of the QIAvac 24. Close unused positions with luer caps and connect QIAvac 24 to a vacuum source.

\* When working with chemicals, always wear a suitable lab coat, disposable gloves, and protective goggles. For more information, please consult the appropriate material safety data sheets (MSDSs) available from the product supplier.

### 3b. QIAvac 6S manifold (see page 31):

Open QIAvac 6S lid. Place QIAvac Luer Adapter(s), or blanks to seal unused slots, into the slots of QIAvac top plate, and close the QIAvac 6S lid. Place the waste tray inside the QIAvac base, and place the top plate squarely over the base. Attach the QIAvac 6S to a vacuum source.

Insert each MinElute column into a luer connector on the Luer Adapter(s) in the manifold. Seal unused luer connectors with plugs provided with the QIAvac Luer Adapter Set.

**Other vacuum manifolds:** follow the suppliers instructions. Insert each MinElute column into a luer connector.

### 4. After the gel slice has dissolved completely, check that the color of mixture is yellow (similar to Buffer QG without dissolved agarose).

**Note:** If the color of the sample is orange or violet, add 10  $\mu$ l of 3 M sodium acetate, pH 5.0, and mix. The color of the mixture will turn to yellow.

The adsorption of DNA to the membrane is efficient only at pH  $\leq 7.5$ . Buffer QG contains a pH indicator which is yellow at pH  $\leq 7.5$  and orange or violet at higher pH, allowing easy determination of the optimal pH for DNA binding.

### 5. Add 1 gel volume of isopropanol to the sample and mix by inverting the tube several times.

For example, if the agarose gel slice is 100 mg, add 100  $\mu$ l isopropanol. Do not centrifuge the sample at this stage.

### 6. To bind DNA, pipet the sample onto the MinElute column and apply vacuum. After the sample has passed through the column, switch off vacuum source.

For maximum recovery, transfer all traces of sample to the column.

The maximum volume of the column reservoir is 800  $\mu$ l. For sample volumes of more than 800  $\mu$ l, simply load again.

### 7. Add 500 $\mu$ l of Buffer QG to the MinElute column and apply vacuum.

### 8. To wash, add 750 $\mu$ l of Buffer PE to the MinElute column and apply vacuum.

**Note:** If the DNA will be used for salt sensitive applications, such as blunt-end ligation and direct sequencing, let the column stand 2–5 min after addition of Buffer PE before applying vacuum.

### 9. Transfer the MinElute column to a clean 1.5 ml microfuge tube or to a provided 2 ml collection tube. Centrifuge for 1 min at $\geq 10,000 \times g$ (~13,000 rpm).

**IMPORTANT:** This spin is necessary to remove residual ethanol (Buffer PE).

### 10. Place the MinElute column in a clean 1.5 ml microcentrifuge tube.

11. To elute DNA, add 10  $\mu$ l of Buffer EB (10 mM Tris-Cl, pH 8.5) or H<sub>2</sub>O to the center of the membrane, let the columns stand for 1 min, and then centrifuge for 1 min.

**IMPORTANT:** Ensure that the elution buffer is dispensed directly onto the center of the membrane for complete elution of bound DNA. The average eluate volume is 9  $\mu$ l from 10  $\mu$ l elution buffer volume.

Elution efficiency is dependent on pH. The maximum elution efficiency is achieved between pH 7.0 and 8.5. When using water, make sure that the pH value is within this range, and store DNA at  $-20^{\circ}\text{C}$  as DNA may degrade in the absence of a buffering agent. The purified DNA can also be eluted in TE\* (10 mM Tris-Cl, 1 mM EDTA, pH 8.0), but the EDTA may inhibit subsequent enzymatic reactions.

\* When working with chemicals, always wear a suitable lab coat, disposable gloves, and protective goggles. For more information, please consult the appropriate material safety data sheets (MSDSs) available from the product supplier.

# MinElute Reaction Cleanup Kit Protocol

## using a microcentrifuge

This protocol is designed to purify double-stranded DNA fragments from all enzymatic reactions e.g., restriction digestion and labeling resulting in high end-concentrations of DNA (see page 10). Fragments ranging from 70 bp to 4 kb are purified from enzymes, primers, nucleotides, and salts using the MinElute Reaction Cleanup columns in a microcentrifuge. The DNA-binding capacity of the MinElute columns is 5 µg.

### Important points before starting

- The yellow color of Buffer ERC indicates a pH  $\leq 7.5$ .
- Add ethanol\* (96-100%) to Buffer PE before use (see bottle label for volume).
- All centrifuge steps are carried out at  $\geq 10,000 \times g$  (~13,000 rpm) in a conventional tabletop microcentrifuge.
- 3 M sodium acetate\*, pH 5.0, may be required.

**1. Add 300 µl of Buffer ERC to the enzymatic reaction and mix. The maximum volume of enzymatic reaction that can be processed per MinElute column is 100 µl.**

If the enzymatic reaction is in a volume of  $< 20 \mu\text{l}$ , adjust the volume to 20 µl. If the enzymatic reaction exceeds 100 µl, split your reaction, add 300 µl of Buffer ERC to each aliquot of the split reaction and use the appropriate number of MinElute columns.

**2. Check that the color of the mixture is yellow (similar to Buffer ERC without the enzymatic reaction).**

If the color of the mixture is orange or violet, add 10 µl of 3 M sodium acetate, pH 5.0, and mix. The color of the mixture will turn to yellow.

**3. Place a MinElute column in a 2 ml collection tube in a suitable rack.**

**4. To bind DNA, apply the sample to the MinElute column and centrifuge for 1 min.**

To obtain maximal recovery, transfer all traces of sample to the spin column.

**5. Discard the flow-through and place the MinElute column back into the same tube.**

**6. To wash, add 750 µl Buffer PE to the MinElute column and centrifuge for 1 min.**

**7. Discard the flow-through and place the MinElute column back in the same tube. Centrifuge the column for an additional 1 min at maximum speed.**

**IMPORTANT:** Residual ethanol from Buffer PE will not be completely removed unless the flow-through is discarded before this additional centrifugation.

\* When working with chemicals, always wear a suitable lab coat, disposable gloves, and protective goggles. For more information, please consult the appropriate material safety data sheets (MSDSs) available from the product supplier.



8. Place the MinElute column in a clean 1.5 ml microcentrifuge tube.
9. To elute DNA, add 10  $\mu$ l Buffer EB (10 mM Tris-Cl, pH 8.5) or H<sub>2</sub>O to the center of the membrane, let the column stand for 1 min, and then centrifuge for 1 min.

**IMPORTANT:** Ensure that the elution buffer is dispensed directly onto the membrane for complete elution of bound DNA. The average volume of eluate is 9  $\mu$ l from 10  $\mu$ l elution buffer. Elution efficiency is dependent on pH. The maximum elution efficiency is achieved between pH 7.0 and 8.5. When using water, make sure that the pH value is within this range, and store DNA at  $-20^{\circ}\text{C}$  as DNA may degrade in the absence of a buffering agent. The purified DNA can also be eluted in TE (10 mM Tris-Cl, 1 mM EDTA, pH 8.0), but the EDTA may inhibit subsequent enzymatic reactions.

\* When working with chemicals, always wear a suitable lab coat, disposable gloves, and protective goggles. For more information, please consult the appropriate material safety data sheets (MSDSs) available from the product supplier.

# MinElute Reaction Cleanup Kit Protocol

## using a vacuum manifold

MinElute columns can also be used on any vacuum manifold with luer connectors, e.g., QIAvac 24 (or QIAvac 6S with Luer Adapters). The following protocol is designed to purify double-stranded DNA from all enzymatic reactions e.g., restriction digestion and labeling resulting in high end-concentrations of DNA (see page 10). Fragments ranging from 70 bp to 4 kb are purified from enzymes, primers, nucleotides, and salts using vacuum-driven sample processing.

### Important points before starting

- The yellow color of Buffer ERC indicates a pH  $\leq 7.5$ .
- Add ethanol\* (96–100%) to Buffer PE before use (see bottle label for volume).
- Switch off vacuum between steps to ensure that a consistent, even vacuum is applied during manipulations.
- 3 M sodium acetate\*, pH 5.0, may be required.

### 1. Add 300 $\mu$ l of Buffer ERC to the enzymatic reaction and mix. The maximum volume of the enzymatic reaction is 100 $\mu$ l.

If the enzymatic reaction is in a volume of  $< 20 \mu$ l, adjust the volume to 20  $\mu$ l. If the enzymatic reaction exceeds 100  $\mu$ l, split the reaction, add 300  $\mu$ l of Buffer ERC to each aliquot of the split reaction and use the appropriate number of MinElute columns.

### 2. Check that the color of the mixture is yellow (similar to Buffer ERC without the enzymatic reaction).

If the color of the mixture is orange or violet, add 10  $\mu$ l of 3 M sodium acetate, pH 5.0, and mix. The color of the mixture will turn to yellow.

### 3. Prepare the vacuum manifold and MinElute columns:

#### 3a. QIAvac 24 (see page 30):

Place the QIAvac 24 lid on top of the QIAvac 24 base. Make sure that the gasket fits tightly in the groove of the QIAvac 24 lid.

Insert up to 24 MinElute columns into the luer extensions of the QIAvac 24. Close unused positions with luer caps and connect QIAvac 24 to a vacuum source.

\* When working with chemicals, always wear a suitable lab coat, disposable gloves, and protective goggles. For more information, please consult the appropriate material safety data sheets (MSDSs) available from the product supplier.

**3b. QIAvac 6S manifold (see page 31):**

Open QIAvac 6S lid. Place QIAvac Luer Adapter(s), or blanks to seal unused slots, into the slots of QIAvac top plate, and close the QIAvac 6S lid. Place the waste tray inside the QIAvac base, and place the top plate squarely over the base. Attach the QIAvac 6S to a vacuum source.

Insert each MinElute column into a luer connector on the Luer Adapter(s) in the manifold. Seal unused luer connectors with plugs provided with the QIAvac Luer Adapter Set.

**3c. Other vacuum manifolds:** follow the supplier's instructions. Insert each MinElute column into a luer connector.

**4. To bind DNA, load the samples into the MinElute columns by decanting or pipetting, and apply vacuum. After the samples have passed through the column, switch off the vacuum source.**

For maximum recovery, transfer all traces of sample to the column.

**5. To wash, add 750  $\mu$ l of Buffer PE to each MinElute column and apply vacuum.**

**6. Transfer each MinElute column to a microcentrifuge tube or the provided 2 ml collection tubes. Centrifuge for 1 min at  $\geq 10,000 \times g$  (~13,000 rpm).**

**IMPORTANT:** This spin is necessary to remove residual ethanol (Buffer PE).

**7. Place each MinElute column into a clean 1.5 ml microcentrifuge tube.**

**8. To elute DNA, add 10  $\mu$ l of Buffer EB (10 mM Tris-Cl, pH 8.5) or H<sub>2</sub>O to the center of each membrane, let the columns stand for 1 min, and then centrifuge.**

**IMPORTANT:** Ensure that the elution buffer is dispensed directly onto the center of the membrane for complete elution of bound DNA. The average eluate volume is 9  $\mu$ l from 10  $\mu$ l elution buffer volume.

Elution efficiency is dependent on pH. The maximum elution efficiency is achieved between pH 7.0 and 8.5. When using water, make sure that the pH value is within this range, and store DNA at  $-20^{\circ}\text{C}$  as DNA may degrade in the absence of a buffering agent. The purified DNA can also be eluted in TE (10 mM Tris-Cl, 1 mM EDTA, pH 8.0), but the EDTA may inhibit subsequent enzymatic reactions.

\* When working with chemicals, always wear a suitable lab coat, disposable gloves, and protective goggles. For more information, please consult the appropriate material safety data sheets (MSDSs) available from the product supplier.

## Troubleshooting Guide

This troubleshooting guide may be helpful in solving any problems which may arise. The scientists at QIAGEN Technical Services are always happy to answer any questions you may have about either the information and protocol(s) in this handbook or molecular biology applications (see back cover).

### Comments and suggestions

---

#### Low or no recovery

- |    |                                      |   |
|----|--------------------------------------|---|
| a) | Buffer PE did not contain ethanol    | Ethanol must be added to Buffer PE (concentrate) before use. Repeat procedure with correctly prepared Buffer PE.  |
| b) | Inappropriate elution buffer         | DNA will only be eluted efficiently in the presence of low-salt buffer (e.g., Buffer EB: 10 mM Tris-Cl, pH 8.5) or water. See "Elution in low-salt solutions", page 13. |
| c) | Elution buffer incorrectly dispensed | Add elution buffer to the center of the MinElute membrane to ensure that the buffer completely covers the membrane.   |

#### Gel

- |    |  |   |
|----|--|---|
| d) | Gel slice incompletely solubilized   | After addition of Buffer QG to the gel slice, mix by vortexing the tube every 2–3 minutes during the 50°C incubation. DNA will remain in any undissolved agarose.   |
| e) | pH of electrophoresis buffer too high (binding mixture turns orange or violet)   | The electrophoresis buffer has been repeatedly used or incorrectly prepared, resulting in a sample pH that exceeds the buffering capacity of Buffer QG and leads to inefficient DNA binding. Add 10 µl of 3 M sodium acetate, pH 5.0 to the sample and mix. The color of the mixture will turn yellow indicating the correct pH for DNA binding. Even for binding mixtures with only small color changes (slight orange color), add the 10 µl sodium acetate. |
| f) | Cloudy and gelatinous appearance of sample mixture after addition of isopropanol | This may be due to salt precipitation, and will disappear upon mixing the sample.   |

**Gel:** refers to MinElute Gel Extraction Kits only

**PCR:** refers to MinElute PCR Purification Kits only

**Reaction Cleanup:** refers to MinElute Reaction Cleanup Kits only

Other notes refer to all kits.

### **Gel/Reaction Cleanup**

- g) Binding mixture turns orange or violet      The pH in the sample exceeds the buffer capacity of Buffer QG or ERC respectively. Add 20  $\mu$ l of 3 M sodium acetate, pH 5.0, to the sample and mix. The color of the mixture will turn yellow indicating the correct pH for DNA binding. Even for samples with slight color changes (orange color), add 10  $\mu$ l sodium acetate.

### **Reaction Cleanup**

- h) Sample volume too high or low      The sample volume must be in the range of 20–100  $\mu$ l.

### **DNA does not perform well, e.g., in ligation reaction**

- a) Salt concentration in eluate too high      Modify the wash step by incubating the MinElute column for 5 min at room temperature after adding 750  $\mu$ l of Buffer PE, then centrifuge.
- b) Eluate contains residual ethanol      Ensure that the wash flow-through is drained from the collection tube and that the MinElute column is then centrifuged at  $\geq 10,000 \times g$  (~13,000 rpm) for 1 min.

### **Gel**

- c) Eluate contaminated with agarose      The gel slice is incompletely solubilized or weighs >400 mg. be sure to vortex the gel slice in Buffer QG every 2–3 minutes during the solubilization step.

### **PCR**


- d) Eluate contains primer-dimers      Primer-dimers are >20 bp, and are not completely removed. After the binding step, wash the MinElute column with 750  $\mu$ l of a 35% (w/v) guanidine hydrochloride solution (35 g in 100 ml). Continue with the Buffer PE wash step and the elution step as in the protocol.
- e) Eluate contains denatured ssDNA, which appears as smaller smeared band on an analytical gel      Use the eluted DNA to prepare the subsequent enzymatic reaction but omit the enzyme. To reanneal the ssDNA, incubate the reaction mixture at 95°C for 2 min, and allow the tube to cool slowly to room temperature. Add the enzyme and proceed as usual. Alternatively, the DNA can be eluted in 10 mM Tris buffer containing 10 mM NaCl. The salt and buffering agent promote the renaturation of DNA strands. However the salt concentration of the eluate must then be considered for subsequent applications.

## Appendix: QIAvac Vacuum Manifolds

### Handling Guidelines for QIAvac 24

The following guidelines should be followed when working with QIAvac 24.

- Always place QIAvac 24 on a secure bench top or work area. If dropped, the QIAvac 24 manifold may crack.
- Always store QIAvac 24 clean and dry. To clean, simply rinse all components with distilled water and allow to air dry or dry with paper towels.
- The components of QIAvac 24 are not resistant to certain solvents (Table 3). If these solvents are spilled on the unit, rinse it thoroughly with water.
- To ensure consistent performance, do not apply silicone or vacuum grease to any part of the QIAvac 24 manifold.
- Always use caution and wear safety glasses when working near a vacuum manifold under pressure.
- Contact QIAGEN Technical Services or your local distributor for information concerning spare or replacement parts.

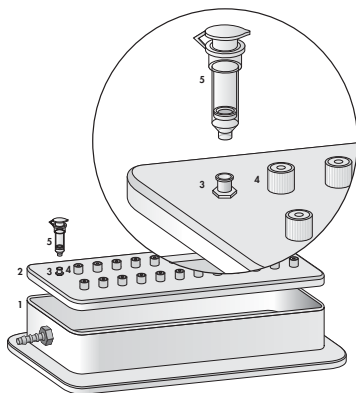

**Figure 5** Components of the QIAvac 24 manifold.

1. QIAvac 24 base
2. QIAvac 24 lid
3. Luer extension of QIAvac 24
4. Luer extension closed with luer cap
5. MinElute column\*

\* Not included with QIAvac 24. Included in appropriate kits.


**Table 3. Chemical Resistance Properties of QIAvac 24**

<b>Resistant to:</b>		<b>Not resistant to:</b>
Chlorine bleach	Acetone	Phenol
Hydrochloric acid	Chromic acid	Benzene
Sodium chloride	Concentrated alcohols	Chloroform
Sodium hydroxide	Chaotropic salts	Ethers
Urea	SDS	Toluene
Acetic acid	Tween 20	

## Handling Guidelines for QIAvac 6S

QIAvac 6S facilitates DNA cleanup with MinElute by providing a convenient modular vacuum manifold, which, in combination with QIAvac Luer Adapters, allows easy processing of MinElute spin columns, as an alternative to centrifugation. The following recommendations should be followed when handling the QIAvac 6S vacuum manifold.

- Always store QIAvac 6S vacuum manifold clean and dry. To clean, simply rinse all components with water and dry with paper towels. Do not air-dry, as the screws may rust and need to be replaced. Do not use abrasives or solvents.
- Always place the QIAvac 6S vacuum manifold on a secure bench top or work area. If dropped, the manifold may crack.
- The components of QIAvac manifolds are not resistant to ethanol, methanol, or other organic solvents (Table 4). Do not bring solvents into contact with the vacuum manifold. If solvents are spilled on the unit, rinse thoroughly with distilled water, and do not incubate acrylic components in alcohol-containing reagents for long periods of time. Ensure that no residual Buffer PE remains in the vacuum manifold.
- To ensure consistent performance, do not apply silicone or vacuum grease to any part of the QIAvac 6S manifold. The spring lock on the top plate and the self-sealing gasket provide an airtight seal when vacuum is applied to the assembled unit. To maximize gasket lifetime, rinse the gasket free of salts and buffers after each use and dry with paper towels before storage.
- Remove blanks from the slots of the top plate after use and store them under the manifold.

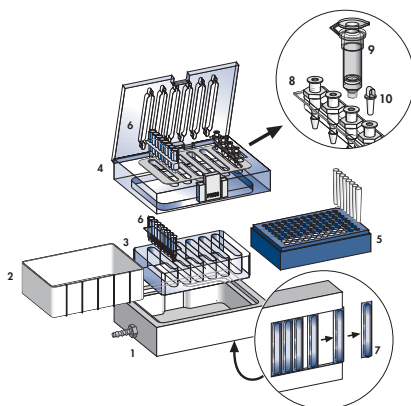

**Figure 6** Components of the QIAvac 6S manifold.

- |   |  |
|---|--|
| 1. QIAvac base, which holds a waste tray, a strip holder, or a microtube rack | 5. Microtube rack                                    |
| 2. Waste tray   | 6. 8-well strip <sup>†</sup>                         |
| 3. QIAvac strip holder to hold 8-well strips                                  | 7. Blanks to seal unused slots                       |
| 4. QIAvac top plate with slots for 8-well strips or QIAvac Luer Adapters      | 8. QIAvac Luer Adapter <sup>‡</sup>                  |
|   | 9. MinElute column <sup>§</sup>                      |
|   | 10. Plug to seal unused luer connectors <sup>†</sup> |

\* MinElute kits are not available in 8-well strip format.

<sup>†</sup> Not included with QIAvac 6S. Must be purchased separately.

<sup>‡</sup> Not included with QIAvac 6S. Included in appropriate kits.

**Table 4. Chemical Resistance Properties of QIAvac 6S**

Resistant to:		Not resistant to:
Chlorine bleach (12%)	Acetic acid	Benzene
Hydrochloric acid	Acetone	Chloroform
Sodium chloride	Chromic acid	Ethers
Sodium hydroxide	Phenol	Toluene
Urea	Concentrated alcohols	

## References

1. Vogelstein, B. and Gillespie, D. (1979) Preparative and analytical purification of DNA from agarose. *Proc. Natl. Acad. Sci. USA* **76**, 615.
2. Hamaguchi, K. and Geiduschek, E.P. (1962) The effect of electrolytes on the stability of deoxyribonucleate helix. *J. Am. Chem. Soc.* **84**, 1329.



# Ordering Information

Product	Contents	Cat. no.
<b>MinElute Kits</b>		
MinElute PCR Purification Kit (50)	50 MinElute Spin Columns, Buffers, Collection Tubes (2 ml)	28004
MinElute PCR Purification Kit (250)	250 MinElute Spin Columns, Buffers, Collection Tubes (2 ml)	28006
MinElute Gel Extraction Kit (50)	50 MinElute Spin Columns, Buffers, Collection Tubes (2 ml)	28604
MinElute Gel Extraction Kit (250)	250 MinElute Spin Columns, Buffers, Collection Tubes (2 ml)	28606
MinElute Reaction Cleanup Kit (50)	50 MinElute Spin Columns, Buffers, Collection Tubes (2 ml)	28204
MinElute Reaction Cleanup Kit (250)	250 MinElute Spin Columns, Buffers, Collection Tubes (2 ml)	28206
<b>Individual Buffers</b>		
Buffer PB (500 ml)	500 ml Binding Buffer	19066
Buffer PE (concentrate, 100 ml)	100 ml Wash Buffer (5x concentrate for 500 ml buffer)	19065
Buffer QG* (250 ml)	250 ml Solubilization and Binding Buffer (with pH indicator)	19063

\* Additional Buffer QG may be required for routine purifications from gel slices >300 mg from gels containing >2% agarose.

## Ordering Information

Product	Contents	Cat. no.
<b>QIAvac Manifolds and Accessories</b>		
QIAvac 24	Vacuum manifold for processing 1–24 spin columns: includes QIAvac 24 Base, Lid, Luer Caps	19403
QIAvac 6S	Vacuum manifold for processing 1–24 spin columns: includes Top Plate with flip-up lid, Base, Waste Tray, Blanks, Strip Holder, Rack of Collection Microtubes (1.2 ml)	19503
QIAvac Luer Adapter Set*	For processing 1–24 QIAGEN spin columns on QIAvac 6S: 6 adapters with 4 luer connectors each, 24 plugs	19541

\* Compatible only with QIAvac Top Plates containing flip-up lid.

# QIAGEN Companies

Please see the back cover for contact information for your local QIAGEN office.

## QIAGEN Distributors

### Argentina

Tecnolab S.A.  
Tel: (011) 4555 0010  
Fax: (011) 4553 3331  
E-mail: info@tecnolab.com.ar  
Web site: www.tecnolab.com.ar

### Austria

VWR International GmbH  
Tel: (01) 576 00 0  
Fax: (01) 576 00 600  
E-mail: info@at.vwr.com  
Web site: www.vwr.com

### Brazil

Uniscience do Brasil  
Tel: 011 3622 2320  
Fax: 011 3622 2323  
E-mail: info@uniscience.com  
Web site: www.uniscience.com

### China

Gene Company Limited  
Tel: (852)2896-6283  
Fax: (852)2515-9371  
E-mail: info@gench.com  
Hong Kong:  
info@gench.com  
Beijing:  
info\_bj@genecompany.com  
Shanghai:  
info\_sh@genecompany.com  
Chengdu:  
gene@public.cd.sc.cn  
Guangzhou:  
info\_gz@genecompany.com

### Croatia

INEL Medicinska Tehnika d.o.o.  
Tel: (01) 2984-898  
Fax: (01) 6520-966  
E-mail: inel-medicinska-tehnika@zg.htnet.hr

### Cyprus

Scientronics Ltd  
Tel: 02-357 22 765416  
Fax: 02-357 22 764614  
E-mail: a.sarpetsas@biotronics.com.cy

### Czech Republic

BIO-CONSULT spol. s.r.o.  
Tel/Fax: (420) 2 417 29 792  
E-mail: info@bioconsult.cz  
Web site: www.bioconsult.cz

### Denmark

VWR International A/S  
Tel: 43 86 87 88  
Fax: 43 86 87 90  
E-mail: info@dk.vwr.com  
Web site: www.vwr.com

### Egypt

Clinilab  
Tel: 52 57 212  
Fax: 52 57 210  
E-mail: Clinilab@link.net

### Finland

VWR International Oy  
Tel: (09) 804 551  
Fax: (09) 8045 5200  
E-mail: info@fi.vwr.com  
Web site: www.vwr.com

### Greece

BioAnalytica S.A.  
Tel: (210)-640 03 18  
Fax: (210)-646 27 48  
E-mail: bioanalyt@hol.gr  
Web site: www.bioanalytica.gr

### Hungary

Kasztel-Med Co. Ltd.  
Tel: (01) 385 3887  
Fax: (01) 381 0695  
E-mail: info@kasztel.hu  
Web site: www.kasztel.hu

### India

Genetix  
Tel: (011)2542 1714  
or (011)2515 9346  
Fax: (011)2546 7637  
E-mail: genetix@nda.vsnl.net.in

### Israel

Westburg (Israel) Ltd.  
Tel: 08-6900655  
or 1-800 20 22 20 (toll free)  
Fax: 08-6900650  
E-mail: info@westburg.co.il  
Web site: www.westburg.co.il

### Korea

LRs Laboratories, Inc.  
Tel: (02) 924-86 97  
Fax: (02) 924-86 96  
E-mail: webmaster@lrslab.co.kr  
Web site: www.lrslab.co.kr

### Malaysia

RESEARCH BIOLABS SDN. BHD.  
Tel: (603)-8070 3101  
Fax: (603)-8070 5101  
E-mail: biolabs@tm.net.my  
Web site: www.researchbiolabs.com

### Mexico

Quimica Valaner S.A. de C.V.  
Tel: (55) 55 25 57 25  
Fax: (55) 55 25 56 25  
E-mail: ventas@valaner.com  
Web site: www.valaner.com

### New Zealand

Biolab Ltd  
Tel: (09) 980 6700  
or 0800 933 966  
Fax: (09) 980 6788  
E-mail: biosciences@nzl.biolabgroup.com  
Web site: www.biolabgroup.com/nzl

### Norway

VWR International AS  
Tel: 02 29 0  
Fax: 22 90 00 40  
E-mail: info@no.vwr.com  
Web site: www.vwr.com

### Poland

Syngen Biotech Sp.z.o.o.  
Tel: (071) 351 41 06  
or 0601 70 60 07  
Fax: (071) 351 04 88  
E-mail: info@syngen.pl  
Web site: www.syngen.pl

### Portugal

IZASA PORTUGAL, LDA  
Tel: (21) 424 7312  
Fax: (21) 417 2674  
E-mail: consulatbiotec@izasa.es

### Singapore

Research Biolabs Pte Ltd  
Tel: 67775366  
Fax: 67785177  
E-mail: sales@researchbiolabs.com  
Web site: www.researchbiolabs.com

### Slovak Republic

BIO-CONSULT Slovakia spol. s.r.o.  
Tel/Fax: (02) 5022 1336  
E-mail: bio-cons@edicon.sk  
Web site: www.bioconsult.cz

### Slovenia

MEDILINE d.o.o.  
Tel: (01) 830-8040  
Fax: (01) 830-8070  
or (01) 830-8063  
E-mail: mediline@sicl.net

### South Africa

Southern Cross Biotechnology (Pty) Ltd  
Tel: (021) 671 5166  
Fax: (021) 671 7734  
E-mail: info@scb.co.za

### Spain

IZASA, S.A.  
Tel: (93) 902.20.30.90  
Fax: (93) 902.22.33.66  
E-mail: consulatbiotec@izasa.es

### Sweden

VWR International AB  
Tel: (08) 621 34 00  
Fax: (08) 760 45 20  
E-mail: info@se.vwr.com  
Web site: www.vwr.com

### Taiwan

TAIGEN Bioscience Corporation  
Tel: (02) 2880 2913  
Fax: (02) 2880 2916  
E-mail: taigen@ms10.hinet.net

### Thailand

Theera Trading Co. Ltd.  
Tel: (02) 412-5672  
Fax: (02) 412-3244  
E-mail: theeatrad@samart.co.th

### Turkey

Medek Medikal Ürünler ve Sağlık Hizmetleri A. S.  
Tel: (216) 302 15 80  
Fax: (216) 302 15 88  
E-mail: makialp@med-ek.com

### All other countries

QIAGEN GmbH, Germany

**Australia** = **QIAGEN Pty Ltd** = PO Box 641 = DONCASTER VIC 3108

Orders 03-9840-9800 = Fax 03-9840-9888 = Technical 1-800-243-066

**Belgium** = **QIAGEN Benelux B.V.** = Spoorstraat 50 = 5911 KJ Venlo = The Netherlands

Orders 0800-79612 = Fax 0800-79611 = Technical 0800-79556

**Canada** = **QIAGEN Inc.** = 2800 Argentia Road = Unit 7 = Mississauga = Ontario = L5N 8L2

Orders 800-572-9613 = Fax 800-713-5951 = Technical 800-DNA-PREP (800-362-7737)

**France** = **QIAGEN S.A.** = 3 avenue du Canada = LP 809 = 91974 COURTABOEUF CEDEX

Orders 01-60-920-920 = Fax 01-60-920-925 = Technical 01-60-920-930

**Germany** = **QIAGEN GmbH** = QIAGEN Strasse 1 = 40724 Hilden

Orders 02103-29-12000 = Fax 02103-29-22000 = Technical 02103-29-12400

**Italy** = **QIAGEN S.p.A.** = Via Grosio, 10/10 = 20151 Milano

Orders 02-33430411 = Fax 02-33430426 = Technical 02-33430414

**Japan** = **QIAGEN K.K.** = Forefront Tower II = 13-1, Kachidoki 3 Chome = Chuo-ku, Tokyo 104-0054

Telephone 03-5547-0811 = Fax 03-5547-0818 = Technical 03-5547-0811

**Luxembourg** = **QIAGEN Benelux B.V.** = Spoorstraat 50 = 5911 KJ Venlo = The Netherlands

Orders 8002-2076 = Fax 8002-2073 = Technical 8002-2067

**The Netherlands** = **QIAGEN Benelux B.V.** = Spoorstraat 50 = 5911 KJ Venlo

Orders 0800-0229592 = Fax 0800-0229593 = Technical 0800-0229602

**Switzerland** = **QIAGEN AG** = Auf dem Wolf 39 = 4052 Basel

Orders 061-319-30-30 = Fax 061-319-30-33 = Technical 061-319-30-31

**UK and Ireland** = **QIAGEN Ltd.** = QIAGEN House = Fleming Way = Crawley = West Sussex, RH10 9NQ

Orders 01293-422-911 = Fax 01293-422-922 = Technical 01293-422-999

**USA** = **QIAGEN Inc.** = 27220 Turnberry Lane = Suite 200 = Valencia = CA 91355

Orders 800-426-8157 = Fax 800-718-2056 = Technical 800-DNA-PREP (800-362-7737)

